




Please see Technical Manual TM295 for detailed safety and operating instructions.

INSTRUCTIONS FOR USE OF PRODUCT AS2000.
Configured in SEV.


1

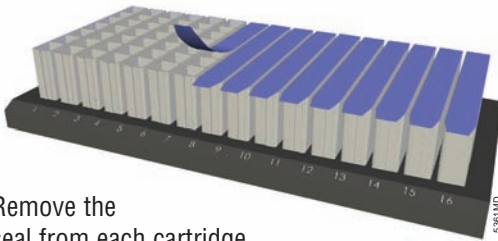

Remove the seal from each cartridge.


5

If you have less than 16 samples,

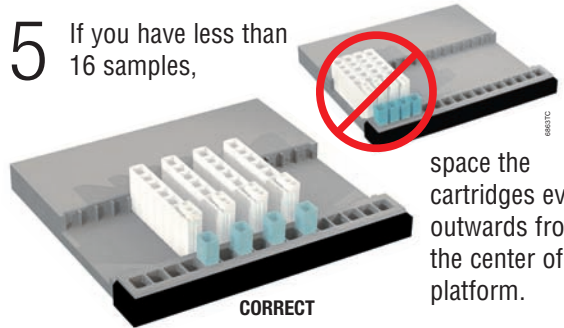

space the cartridges evenly outwards from the center of the platform.

2

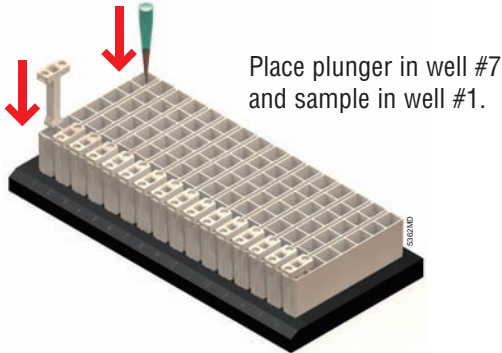

Place plunger in well #7 and sample in well #1.

6




Close the door. Press "Run" to perform purification.

3



Open the door. Press "Run/Stop" to extend platform.

7

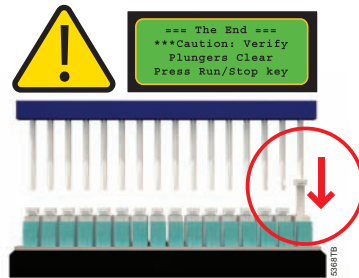


Run completed. Verify that plungers are clear of magnetic rod assembly.

4



Transfer cartridges to platform. Add elution tubes and buffer.

8

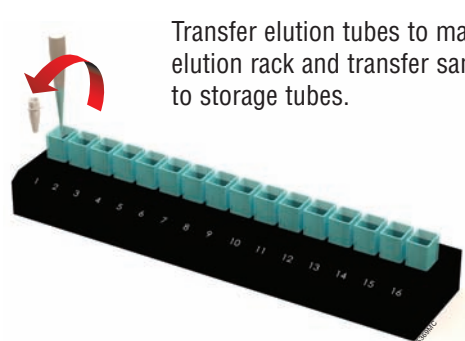

Transfer elution tubes to magnetic elution rack and transfer samples to storage tubes.

ORDERING/TECHNICAL INFORMATION:

www.promega.com • Phone 608-274-4330 or 800-356-9526 • Fax 608-277-2601

©2007 Promega Corporation. All Rights Reserved.

Printed in USA 8/07. 15504-PM-IN. Part# FB086



Promega